

# Uniblitz® CS25

## 25mm Uni-Stable Optical Shutter

### Overview

The Uniblitz CS25 has been designed to provide accurate, repeatable exposures for a wide variety of applications including microscopy, video imaging, and more. The slim form-factor provides a 25mm aperture that can be inserted into a 2.37 inch diameter housing. The CS25 is available in a housed or an un-housed configuration for OEM applications (or simply where spatial limitations are a consideration).

Uni-stable shutter devices, like the CS25, require power to hold the blades in the open state.

### Key Features

- 25mm aperture
- 510A cable included
- Configured for the [VCM-D1 Shutter Driver](#)
- **RoHS Compliant**
- Transfer time on opening:  
9.0 milliseconds
- Transfer time on closing:  
12.0 milliseconds

### Specifications

#### Electrical Specifications

Coil resistance	12 OHMS
Voltage to Open	+65 VDC
Hold Voltage (Nominal) <sup>1</sup>	+5 VDC

#### Mechanical Specifications

Weight Unhoused	36.3 g
Weight Housed	89.6 g
Operating Temp.	0 - 80 °C
Max. Opening Bounce	15%
Max. Closing Bounce	5%
Max. Freq. of Operation <sup>2</sup>	5 Hz / 30 Hz
Number of Shutter Blades	5

<sup>1</sup> Voltage level required across actuator coil when being held in the open position.

<sup>2</sup> (Continuous/Burst) Continuous frequency rating specified at shutter's minimum exposure pulse. Burst frequency rating specified for four (4)




**Need Support?** Please visit our website or email us at [info@uniblitz.com](mailto:info@uniblitz.com).

Tel: +1 585-385-5930 | Fax: 585-385-6004 | 803 Linden Ave. Rochester, NY 14625

Updated 11/21 | Document Version 6.1 | ©2021 Vincent Associates

## Shutter Timing Data

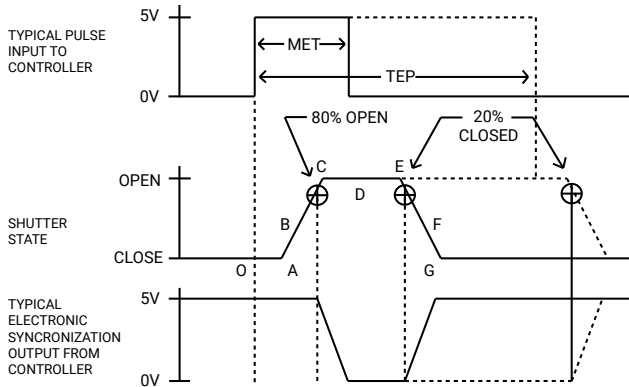

<sup>1</sup> Under no circumstances should any type of lubricant be applied to the shutter blade area. Lubricating the shutter blades will likely slow the shutter down and may eventually render it inoperable.

### CS25 (w/ VCM-D1 and "T" blades) <sup>1</sup>

### Time (msec.)

Parameter	Description	Time (msec.)
O - A	Delay time on opening after current applied	3.0
A - C	Transfer time on opening	9.0
O - C	Total opening time	12.0
C - E	Min. dwell time with min. input pulse	6.0
B - F	Min. equivalent exp. time	17.5
E - G	Transfer time on closing	14.0
A - G	Total window time	29.0
MET	Min. exposure time	15.0
TEP	Typical exposure pulse	>15.0

## Product Options

CS25 **2** **3** **4** **5** **6** - **7** - **8**

Ex: CS25S3T0-EC-21

### 1 Shutter Series:

CS25

### 2 Voltage:

S: Use with VCM-D1

E: Use with D880C or VED24

### 3 Housing:

1: Un-Housed

3: #3 Housing

### 4 Blades: <sup>2</sup>

T: Low Energy (Teflon®)

ZM: High Energy (AlMgF<sub>2</sub>) <sup>3</sup>

### 5 Electronic Sync:

0: Omitted

1: Included

### 6 Connector:

L: 18" Flying Leads

Leave blank for 5-pin Switchcraft connector

### 7 Encapsulated Coil:

EC: Included

Leave blank if not required

### 8 Mount: <sup>4</sup>

21: Zeiss Axiovert    105: C-Mount (Male)

24: Olympus BX/IX    106: C-Mount (Female)

27: Nikon    Leave blank if not required


<sup>2</sup> Other blade coating options may be available by special order.

<sup>3</sup> Input side only; Teflon® coating is on opposite side to protect shutter blade surface. Light source must be input to the reflective side only.

<sup>4</sup> Mounts are only compatible if #3 housing is optioned as well.


## Electronic Sync.

The electronic synchronization system provides a feedback signal (through the driver utilized) after the shutter transfers to the open state. The system incorporates an infrared emitting diode, an infrared sensitive detecting transistor, and an interrupting vane. The vane is attached to the shutter so as to block the light path between the emitter and detector in the closed position. When the shutter transfers to the 80% open position, the vane is removed from the infrared light path, allowing the emitter to switch the detector to the active state. For the CS25, this system uses a similar activation flag attached to the mechanism, which triggers a reflective emitter/detector device. No connection to the designated synchronization pins when no electronic sync. is selected.

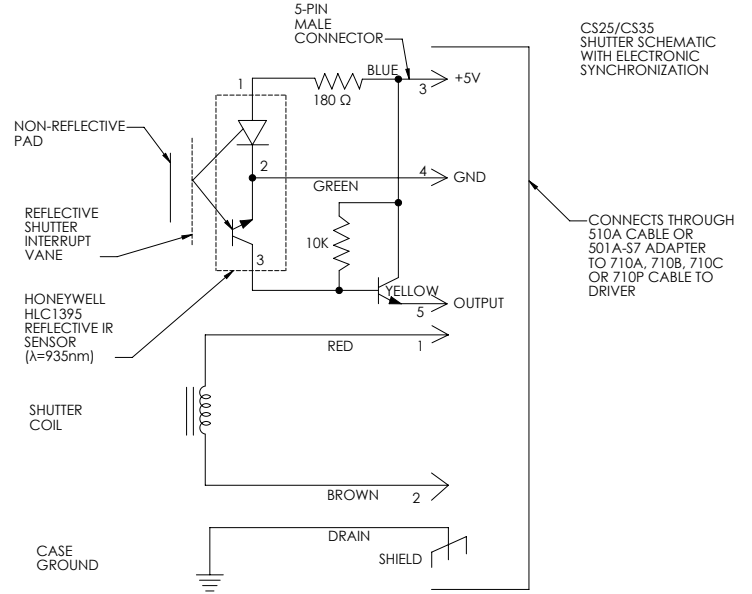


# Uniblitz® CS25 Technical Drawings

## Un-Housed

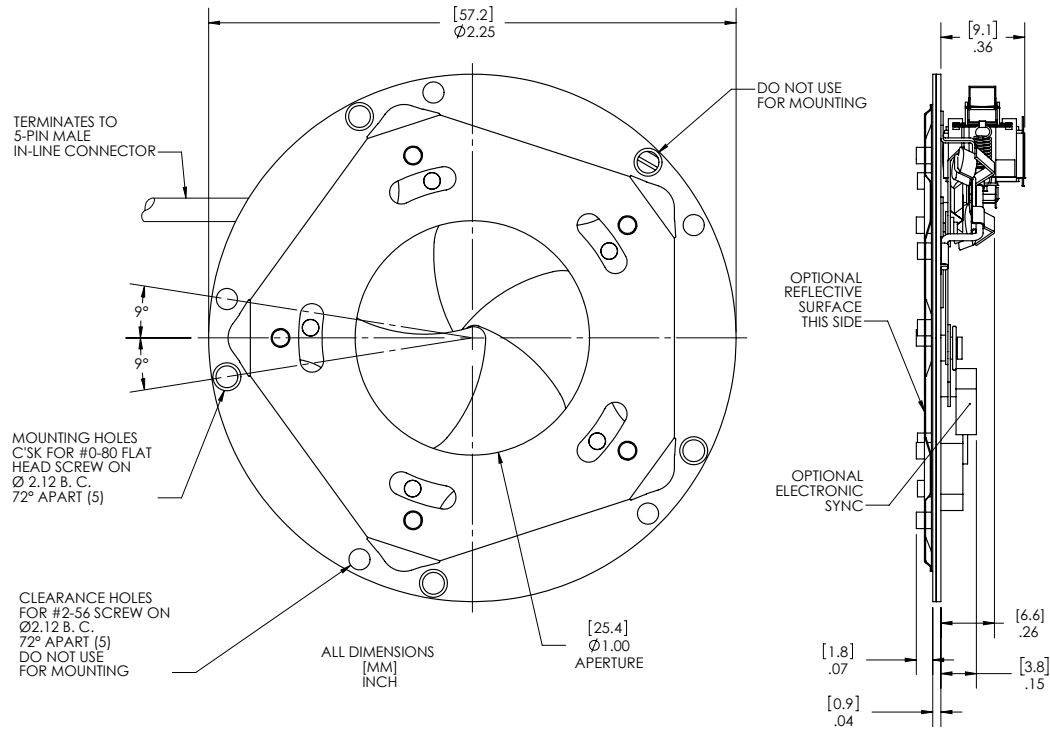


## Housed

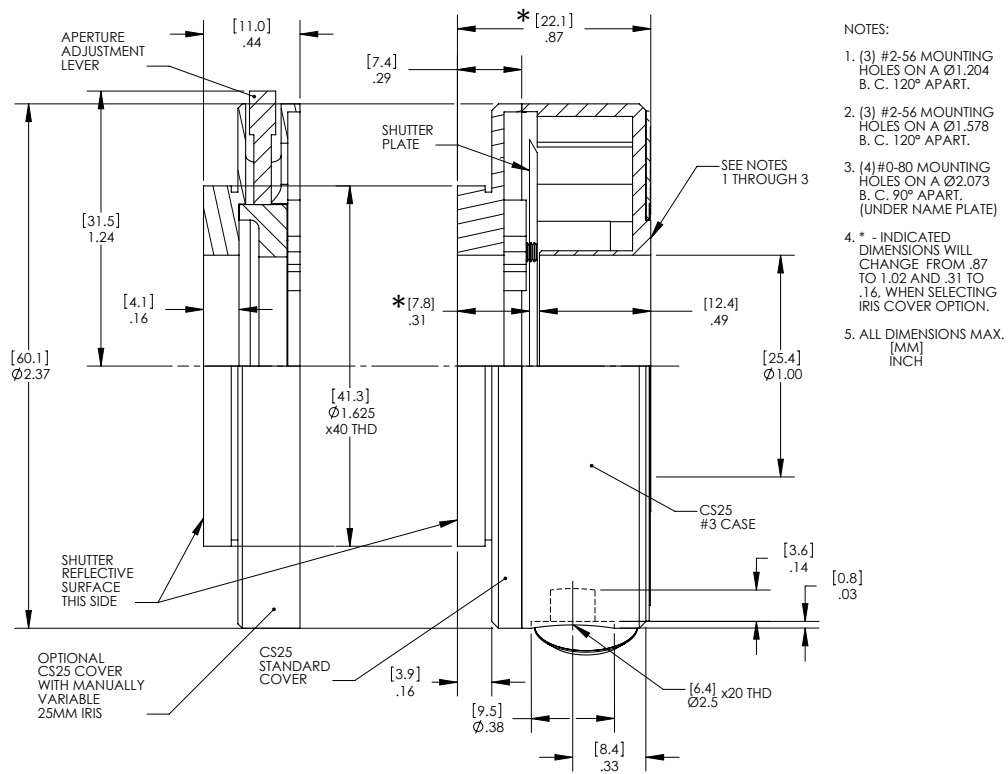

# Uniblitz® CS25 Technical Drawings

## Connector Layout

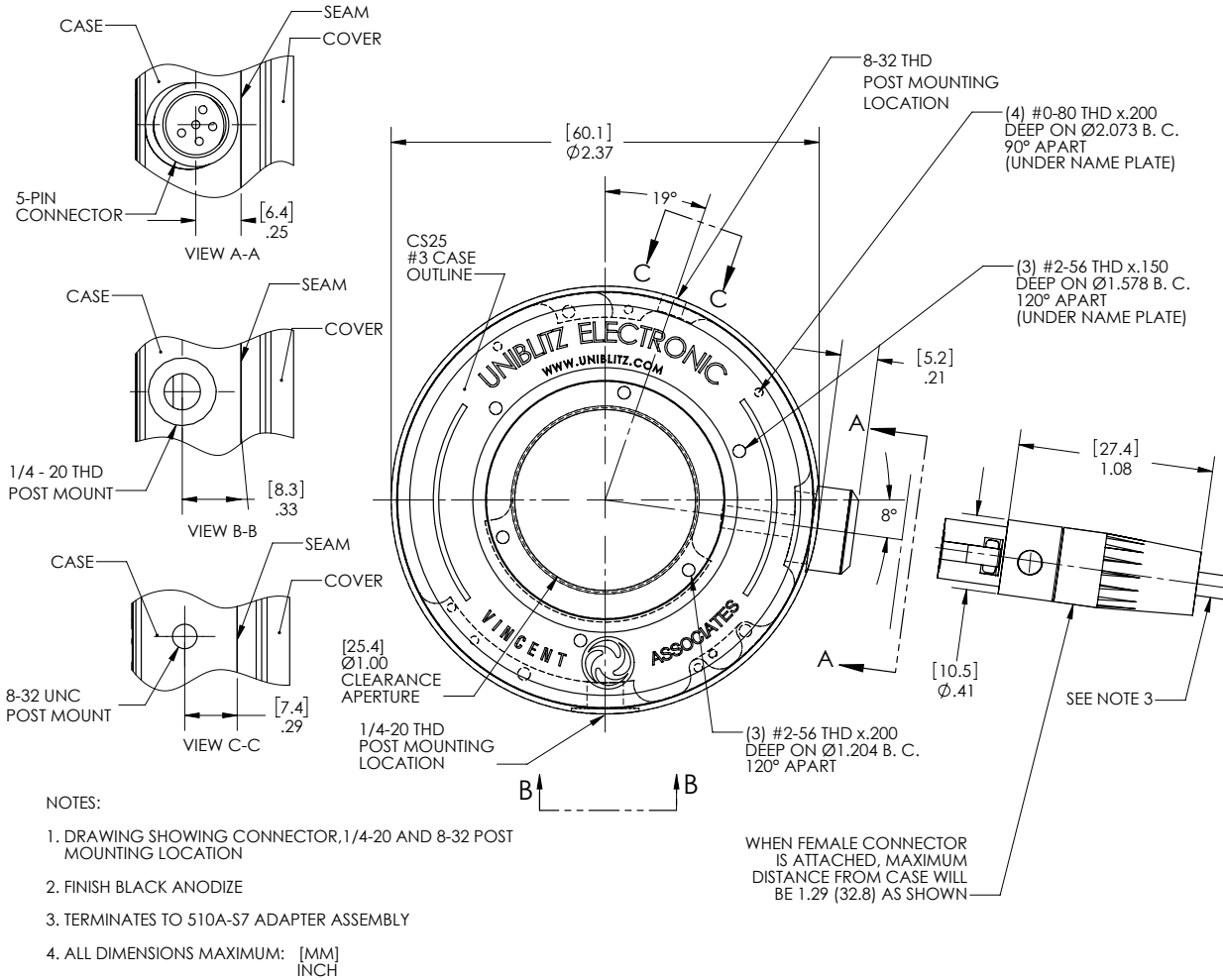

## 510A Interconnect Cable (Included)

