Uniblitz® FS35

35mm Uni-Stable Optical Shutter

Overview

The Uniblitz FS series shutters are designed and optimized to operate directly from +5VDC and do not require a separate driver. Removing the +5VDC (0VDC) closes these shutters. A simple control circuit can be used to operate these shutter devices from a TTL trigger pulse. This control can be also accomplished with our new **VLM1 TTL control interface module**, which is now available.

This low-cost innovation provides the reliability of Uniblitz shutters (typical lifetime >300K cycles) at a single operating voltage.

Key Features

- 35mm aperture
- Default closed operation
- RoHS Compliant
- Transfer time on opening:
 20.0 milliseconds
- Transfer time on closing:
 18.0 milliseconds

Specifications


Electrical Specifications	
Coil resistance	12 OHMS
Voltage to Open	+5 VDC at 425 mA
Hold Voltage	+5 VDC

Mechanical Specifications	
Weight Unhoused	48.5 g
Weight Half-Housed	116.0 g
Weight Housed	183.5 g
Operating Temp.	-40 - +65 C
Max. Opening Bounce	15%
Max. Closing Bounce	5%
Max. Freq. of Operation ¹	12 Hz
Number of Shutter Blades	5

¹ (Continuous/Burst) Continuous frequency rating specified at shutter's minimum exposure time (MET).



Shutter Timing Data

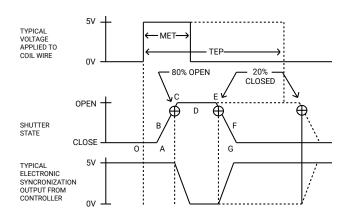

¹ Under no circumstances should any type of lubricant be applied to the shutter blade area. Lubricating the shutter blades will likely slow the shutter down and may eventually render it inoperable.

FS35 (w/ 5VDC and C-PET blades) ¹ Time (msec.)		msec.)
O - A	Delay time on opening after current applied	12.0
A - C	Transfer time on opening	20.0
O - C	Total opening time	32.0
C - E	Min. dwell time with min. input pulse	14.0
B - F	Min. eqivalent exp. time	33.0
E - G	Transfer time on closing	18.0
A - G	Total window time	52.0
MET	Min. exposure time	40.0
TEP	Typical exposure pulse	>40.0

Product Options

FS35 **23456** - **7**

3 Housing: 2 4 Blades: 3 1 Shutter Series: 2 Voltage: FS35: Normally Closed S: Standard 5 VDC 1: Un-Housed **C:** Carbon impregnated plastic (C-PET) FSR35: Normally Open 2: Half-Housed T: Low Energy (Teflon®) 3: Fully-Housed ZM: High Energy (AIMgF2) 4 **6** Connector: **5** Electronic Sync: **7** Encapsulated Coil: 0: Omitted L: 18" Flying Leads EC: Included 5 1: Included Leave blank if not required

Ex: FS35S2C0L-EC

² Various mounting methods available depending on housing option.


³ Other blade coating options may be available by special order.

⁴ Input side only; Teflon® coating is on opposite side to protect shutter blade surface. Light source must be input to the reflective side only.

⁵ Vacuum compatibility up to 10E-10 Torr.

Electronic Sync.

The synchronization system for FS shutter devices incorporates a small magnet mounted to the driving mechanism and a Hall effect sensor. When the device achieves approximately 80% of full open, the magnet causes the Hall effect sensor to change state, producing a signal to indicate that the shutter has switched to the active state. Shown to the left is the FS series shutter schematic which incorporates the electronic synchronization system. There is no connection to the designated synchronization pins when an electronic sync. is not selected.

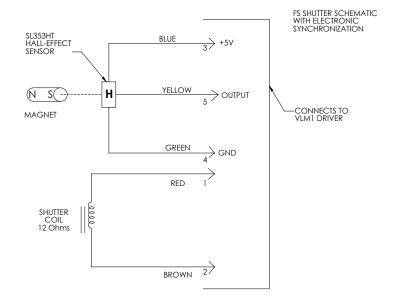

Typical Control Circuit


This simple control circuit to the right can be used to operate the shutter device from a TTL trigger pulse.¹ This control can also be accomplished with the <u>VLM1 interface</u> module.

¹ Note that the shutter will not operate directly from a TTL signal.

+5V. 0.5A 1. ALL RESISTORS 1/8W, 5%, EXCEPT AS NOTED. Red Wire 2. DIODES D1 & D2 PREVENT INPUT VOLTAGES EXCEEDING COIL (A) +5V AND DAMAGING MOSFET Q1. 3. +12V MAXIMUM INPUT FOR 1/8W RESISTORS. SHUTTER COIL DCR = 12 +5V Brown Wire COIL (B) D1 1N4148W 太 R1 Q1 TTL INPUT DMN2046U 1K OR EQUIVALENT D2 R2 本 1N4148W 10K

Uniblitz® FS35 Technical Drawings

Half-Housing - FS35S2C0L

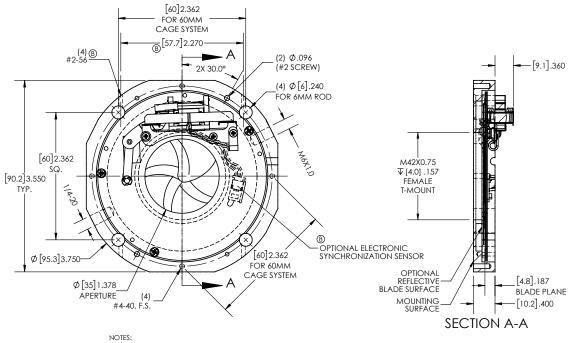


UNITS: [MM] INCH.

LIMITED HIDDEN LINES SHOWN FOR CLARITY.

LEADS WIRES REMOVED FOR CLARITY.

Un-Housed - FS35S1C0L

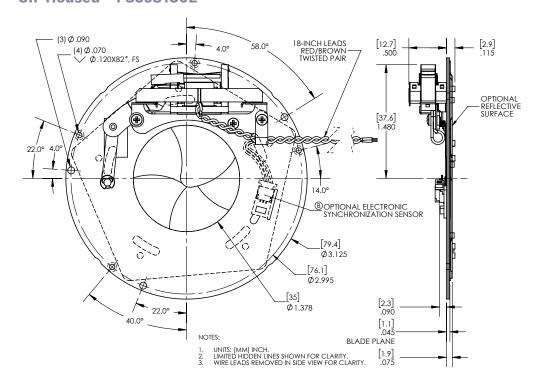

Uniblitz® FS35 Technical Drawings

Full-Housing - FS35S3C0L

